

ANALYSIS OF FIXED GASES FROM UNDERGROUND WATER

Technique : Micro-GC

Column : Tamis

Carrier gas : Helium


Injector : Variable volume

Column temperature : 35°C

Injection time : 300 ms

Column pressure : 35 psi

μTCD sensitivity : High

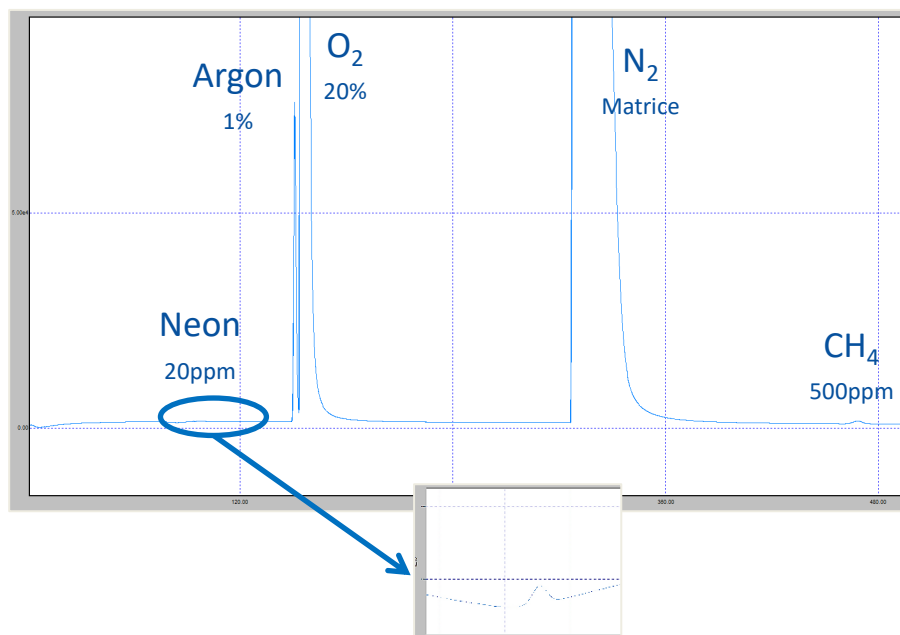

Technique : Micro-GC

Column : PoraPlot Q 8m

Carrier gas : Helium


Injector : Variable volume

Column temperature : 35°C

Injection time : 150 ms

Column pressure : 28 psi

μTCD sensitivity : High

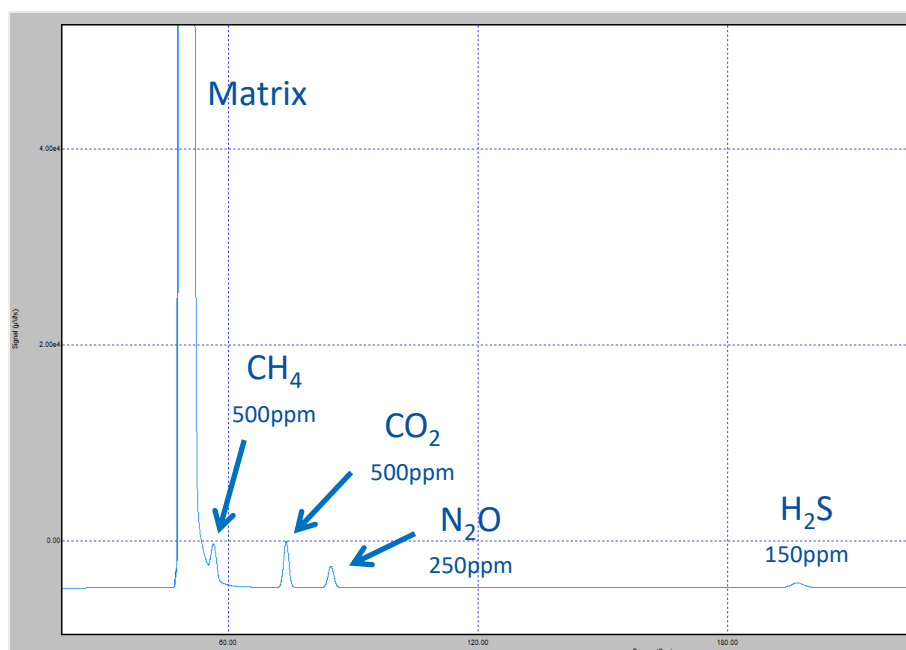

More information : www.sra-instruments.com